ISE F- 9655C EN

Refer to the meter manual for more information.
Caution
 ⚠ Personal injury hazard. Glass components can break. Handle with care to prevent cuts.
 ⚠ Chemical exposure hazard. Refer to the current material safety data sheets (MSDS) for safety protocols.

Introduction
 This probe is a combination fluoride selective electrode with a non-refillable reference. The probe measures fluoride concentration in aqueous solutions.

Technical information

Range	0.01 mg/L (1 x 10 ⁻⁶) to 19,000 mg/L (1 M) F ⁻
pH range	4 to 8
Linear range	0.1 mg/L to 19,000 mg/L Fluoride
Slope	59 mV/pF, 90 to 110% at 25 °C (77 °F)
Ionic strength adjuster (ISA)	Hach ISA 2589999 pillow
	Hach TISAB solution 2829017
Operating temperature	5 to 50 °C (41 to 122 °F)
Interferences	Fe ³⁺ , Al ³⁺ ; remove with ISA or TISAB

Preparation

- 1 Connect the probe.**
- 2 Calibration**
Add 1 ISA pillow or 5.0 mL of TISAB per 25 mL of standard.
- 3 Measurement**
Add 1 ISA pillow or 5.0 mL of TISAB per 25 mL of sample.

Maintenance and storage

Cleaning: rinse with distilled water between each standard and each sample.
 Storage: clean and dry the probe and install the sensor protection cap for storage.

Troubleshooting

The ionic strength and temperature of samples and standards must be the same. Make sure that the stir speed is the same for all standards and samples.
 Damage or deposits on the membrane change the response of the electrode. Remove deposits from the membrane and replace the sensor if damaged.

Warranty

6 months for manufacturing faults only. The warranty does not cover normal wear or improper use.

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